

VAN LONDON

Est. 1961



pHoenix Co.



Fluoroborate Ion Selective Electrode



www.VL-PC.com
info@VL-PC.com

Toll Free (800) 522-7920
Local (713) 772-6641



GENERAL INSTRUCTIONS

Introduction

The Van London-pHoenix Company Fluoroborate Ion Selective Electrode is used to quickly, simply, accurately, and economically measure Fluoroborate in aqueous solutions.

Required Equipment

1. An ion meter
2. Fluoroborate Ion Sensing Electrode
3. Plastic syringe, pipets, volumetric flasks, 150 ml beakers

Required Solutions

1. Deionized or distilled water for solution preparation.
2. Ionic Strength Adjuster (ISA), 2M $(\text{NH}_4)_2\text{SO}_4$, Cat# BF4IS01.
3. Van London-pHoenix Company Fluoroborate Standard, 1000 ppm BF_4^- , Cat. No. BF4AS02.
4. Fluoroborate Electrode Reference Fill Solution, Cat# R001044.

GENERAL PREPARATION

Electrode Preparation

Remove the rubber cap covering the electrode tip and the rubber insert covering the filling hole of the reference electrode. Fill the combination electrode with the filling solution shipped with the electrode to a level just below the fill hole. No preparation is required with a sealed reference electrode. Gently shake the electrode downward in the same manner as a clinical thermometer to remove any air bubbles which may be trapped behind the Fluoroborate membrane. Connect the electrode to the proper terminals of the meter as recommended by the meter manufacturer. Store the black shipping cap for later use.

Electrode Slope Check (for Ion meters which display mV)

1. To a clean, dry, 150 ml beaker, add 100 ml of distilled water and 2 ml of ISA. After assuring that the meter is in the millivolt mode, lower the electrode tip into the solution. Stir moderately. Remove air bubbles on the clear plastic membrane by redipping probe.
2. Using a pipet, add 1 ml of 1000 ppm Fluoroborate standard into the solution. Stir moderately. After 1 minute, record the mV reading.
3. Using a pipet, add 10 ml of the 1000 ppm Fluoroborate standard to the beaker. Stir moderately. After 1 minute, record the mV reading.
4. Determine the difference between the two readings. The electrode is operating correctly if a slope difference of 52-59 mV is found, assuming the solution temperature is 25°C. Slope is defined as the change in mV observed when the concentration changes by a factor of 10.

Measurement using an Ion Meter (in the Concentration Mode)

1. By serial dilution of the 1000 ppm Fluoroborate standard, prepare two ammonia standards whose concentration is near the expected sample concentration. (e.g. 10 ppm and 100 ppm) For example, to make a 100 ppm standard, pipet 10 ml of the 1000 ppm standard into a 100 ml volumetric flask and dilute to volume with deionized water. Next to make a 10 ppm standard, pipet 10 ml of the newly-made 100 ppm standard into a 100 ml volumetric flask and dilute to volume with deionized water. A 1 ppm standard is made by further dilution of the 10 ppm standard. Measure out 100 ml of each standard into

individual 150 ml beakers.

2. Assure that the meter is in the concentration mode and set for a 2-point calibration.
3. Lower the electrode tip into the least concentrated solution. Begin stirring at a constant rate. Add 2 ml of ISA to the solution and continue stirring.
4. After 1 minute, adjust the meter to the concentration of the more dilute Fluoroborate standard and fix the value in the memory according to the meter manufacturer's instructions.
5. Rinse the electrode tip with distilled water and blot dry.
6. Lower the electrode tip into the more concentrated solution. Begin stirring at a constant rate. Add 2 ml of ISA to the solution and continue stirring.
7. After 1 minute, adjust the meter to the concentration of the more concentrated Fluoroborate standard and fix the value in the memory according to the meter manufacturer's instructions.
8. Add 100 ml of the sample and 2 ml of ISA in a 150 ml beaker. Lower the electrode tip into the solution. Begin stirring at a constant rate. Ensure that the meter is in sample mode.
9. After 1 minute, read the concentration directly from the meter display.
10. The electrode should be re-calibrated every 2-3 hours. Simply repeat Steps 2-7 above.

Measuring Hints

As Fluoroborate electrodes are used or stored for long periods, they will experience some deterioration in performance and slope errors will increase. By using the meter's calibration controls this error can be corrected. If an electrode is able to be calibrated and is stable and responsive, it is still a functional electrode and may be used in service even though it no longer meets "new" electrode specifications.

All samples and standards should be at the same temperature for precise measurement. A difference of 1°C in temperature will result in approximately a 2% error.

Always rinse the electrode with distilled water and blot dry between measurements. Use a clean, dry tissue to prevent cross-contamination.

Constant, but not violent, stirring is necessary for accurate measurement.

Check the electrode for air bubbles adhering to the membrane surface after immersion in solution. Agitate the electrode gently to remove the air bubbles.

The sensing membrane is normally subject to water uptake and might appear milky. This has no effect on performance.

A slow responding electrode may be caused by interferences to the electrode. To restore proper performance, soak the electrode in distilled water for about 5 minutes to clean the membrane, rinse, and soak in standard solution for about 5 minutes.

All samples must be aqueous and not contain organics which can dissolve in the membrane or extract out the liquid ion exchanger.

Interferences should be absent. If they are present, use the procedures found in the **Interferences** section to remove them.

The pH range for the Fluoroborate ion electrode is 2.5-11. Neutralize samples outside this range with acid or base to bring them in range.

Fluoroborate sample with either high or low pH must be analyzed immediately after preparation or hydrolysis of fluoroborate to BF_3OH^- , $\text{BF}_2(\text{OH})_2^-$, and $\text{BF}(\text{OH})_3^-$ will occur. Convert boric acid or borate ion to fluoroborate by addition of HF.

ELECTRODE CHARACTERISTICS

Reproducibility

Electrode measurements reproducible to $\pm 2\%$ can be obtained if the electrode is calibrated every hour. Factors such as temperature fluctuation, drift, and noise limit reproducibility.

Interferences

Certain anions are electrode interferences and will cause electrode malfunction, drift or measurement errors if present in high enough levels. The level of interfering common anions that will cause a 10% error at three levels of fluoroborate is given in Table 3.

TABLE 3: Concentration of Possible Interferences Causing a 10% Error at Various Levels of Fluoroborate; Background Ionic Strength of 0.12M $(\text{NH}_4)_2\text{SO}_4$.

Interferences

<u>(ppm)</u>	<u>100 ppm</u>	<u>10 ppm</u>	<u>1 ppm</u>
PO_4^{-3}	7.0×10^4	7.0×10^3	7.0×10^2
HPO_4^{-2}	7.1×10^4	7.1×10^3	7.1×10^2
$\text{H}_2\text{PO}_4^{-1}$	7.2×10^4	7.2×10^3	7.2×10^2
Cl^-	1.6×10^4	1.6×10^3	1.6×10^2
CO_3^{-2}	1.7×10^4	1.7×10^3	1.7×10^2
HCO_3^{-1}	1.7×10^4	1.7×10^3	1.7×10^2
NO_3^{-1}	2.9×10^3	2.9×10^2	2.9×10^1
NO_2^{-1}	4.2×10^2	4.2×10^1	4.2
Br^-	7.4×10^2	7.4×10^1	7.4
CN^-	1.2×10^2	1.2×10^1	1.2
ClO_3^{-1}	4.0×10^1	4.0	4.0×10^{-1}
I^-	6.0	6.0×10^{-1}	6.0×10^{-2}
ClO_4^{-1}	5.0×10^{-1}	5.0×10^{-2}	5.0×10^{-3}

Interferences such as phosphate, bromide, iodide and cyanide can be removed by precipitation with 0.5 grams of silver sulfate added to 100 ml of sample. Interferences such as carbonate and bicarbonate can be removed by conversion to carbon dioxide through acidification of the sample to pH 4.5 with sulfuric acid. Nitrite interference can be removed by adding 0.3 grams of sulfamic acid to 100 ml of sample.

The above interference removal procedures require similar treatment of standards as well as samples.

If the electrode is exposed to high levels of interfering ions which cannot be removed, the electrode reading may drift and the response may become sluggish. Restore performance by soaking in distilled water for 30 minutes followed by soaking in fluoroborate standard for 30 minutes.

Temperature Influences

The electrode response will shift and change slope with change in temperature. Standards and samples should be at

the same temperature. A 2% error results with a 1°C temperature change for a 10 ppm solution.

The electrode can be used at temperatures from 0° - 40°C. Room temperature measurements are recommended, since measurements at temperatures quite different from room temperature may require equilibrium times up to one hour.

Electrode Response

Plotting the mV potential against the Fluoroborate concentration results in a straight line with a slope of about 56 mV between 10 ppm and 100 ppm at 25°C. For Fluoroborate concentrations above 10 ppm BF_4^{-1} , the electrode exhibits good time response (95% of total mV reading in one minute or less). Response times are longer below this value.

Limits of Detection

Fluoroborate concentration down to 0.1 ppm Fluoroborate can be measured in neutral solutions. Since sample contamination can be a factor in low level Fluoroborate measurements, care must be taken in making determinations below 1.0 ppm. The upper limit of detection in pure sodium Fluoroborate solutions is 10,000 ppm.

pH Effects

The operating range of the Fluoroborate electrode is from pH 2.5 to pH 11.

Electrode Life

The Fluoroborate electrode will last six months in normal laboratory use. On-line measurement might shorten operational lifetime to several months. In time, the response time will increase and the calibration slope will decrease to the point calibration is difficult and electrode replacement is required.

Since Fluoroborate electrodes have a limited shelf life, it is important to have a backup electrode which is in working condition when required.

Electrode Storage

The Fluoroborate electrode may be stored for short periods of time in 10 ppm Fluoroborate solution. For longer storage (longer than two weeks), rinse and dry the sensing tip and cover the membrane tip with any protective cap shipped with the electrode. The reference portion of the combination electrode (or the outer chamber of the reference electrode) should be drained of filling solution, if refillable, and the rubber insert placed over the filling hole.

TROUBLESHOOTING HINTS

<u>Symptom</u>	<u>Possible Causes</u>	<u>Next Step</u>
Out of Range Reading	defective meter	check meter with shorting strap (see meter instruction manual)
	defective electrode	check electrode operation
	electrodes not plugged in properly	unplug electrodes and reseal

	reference electrode not filled	be sure reference electrode is filled
	air bubble on membrane	remove bubble by re-dipping electrode
	electrodes not in solution	put electrodes in solution
Noisy or Unstable Readings (readings continuously or rapidly changing)	defective meter	check meter with shorting strap
	air bubble on membrane	remove bubble by re-dipping electrode
	defective electrode	replace electrode
	ISA not used	use recommended ISA
	meter or stirrer not grounded	ground meter or stirrer
	electrode exposed to interferences	soak electrode in Fluoroborate standard
	outer filling solution level too low	fill electrode to level just below the fill hole
Drift (reading slowly changing one direction)	samples and standards at different temperatures	allow solutions to come to room temperature before measurement
	electrode exposed to interferences	soak electrode in Fluoroborate standard
	incorrect reference filling solution	use recommended filling solution
Low Slope or No Slope	standards contaminated or incorrectly made	prepare fresh standards
	ISA not used	use recommended ISA
	standard used as ISA	use ISA
	electrode exposed to interferences	soak electrode in Fluoroborate standard
	defective electrode	check electrode operation
	air bubble on membrane	remove bubble by re-dipping probe
	Incorrect Answer	incorrect standards

wrong units used

apply correct conversion factor:
 $10^{-3}\text{M} = 86.8 \text{ ppm as } \text{BF}_4^{-1}$

sample carryover

rinse electrodes thoroughly between samples

SPECIFICATIONS

Concentration Range:	10,800 ppm to 0.5 ppm
pH Range:	2.5 to 11
Slope:	52 – 59 mV between 10 ppm and 100 ppm at 25°C
Temperature Range:	0° to 40°C
Interferences:	ClO_4^{-1} , I^{-1} , CN^{-1}
Reproducibility:	$\pm 2\%$
Size:	110 mm length 12 mm diameter 1 m cable length
Storage:	store in dilute Fluoroborate standard
Samples:	aqueous solutions only no organic solvents

ELECTRODE THEORY

Electrode Operation

The Fluoroborate electrode consists of an electrode body containing a liquid internal filling solution in contact with a gelled organophilic membrane containing a Fluoroborate ion exchanger. When the membrane is in contact with a solution containing free Fluoroborate ions, an electrode potential develops across the membrane. This electrode potential is measured against a constant reference potential, using a standard pH/mV meter or an ion meter. The level of Fluoroborate ions, corresponding to the measured potential, is described by the Nernst equation:

$$E = E' - S \log X$$

where: E = measured electrode potential
E' = reference potential (a constant)
S = electrode slope (56 mV/decade)
X = level of Fluoroborate ions in solution

The activity, X, represents the effective concentration of the ions in solution. The total Fluoroborate ion concentration, C_t , is the sum of free Fluoroborate ion, C_f , and complexed or bound Fluoroborate ion, C_b . The electrode is able to respond to only the free ions, whose concentration is :

$$C_f = C_t - C_b$$

Since Fluoroborate ions form very few stable complexes, the free ion concentration may be equated to the total ion concentration.

The activity is related to the free ion concentration, C_f , by the activity coefficient, γ , by:

$$X = \gamma C_f$$

Activity coefficients vary, depending on total ions strength, I , defined as:

$$I = 1/2 \sum C_x Z_x^2$$

where: C_x = concentration of ion X
 Z_x = charge of ion X
 \sum = sum of all of the types of ions in the solution

In the case of high and constant ionic strength relative to the sensed ion concentration, the activity coefficient, γ , is constant and the activity, X , is directly proportional to the concentration.